



CARBON / HYDROGEN / SULFUR ANALYZER

CHS-580A

**ELTRA's CHS-580A is the ideal elemental analyzer for simultaneous determination of carbon, hydrogen and sulfur in organic samples. It features a resistance furnace with vertical ceramic tube and uses ceramic crucibles as sample carriers.**

An automatic sample loader allows for analysis of large numbers of samples without operator intervention. The sample loader is available in two versions to accommodate 36 or 130 crucibles. For reliable analysis of acidified samples (TOC analysis), both versions are available in an acid-resistance configuration.

The temperature can be set in steps of 1 °C to a maximum of 1,550 °C. Thanks to sample weights of up to 500 mg and more even inhomogeneous materials can be reliably analyzed. The analyzer can be equipped with up to four independent infrared cells for measuring carbon and sulfur concentrations, allowing for a great variety of applications.

## **BENEFITS CHS-580A**

- | Fast & simultaneous elemental analysis of C, H, S, TOC and TIC
- | Configurable for single element (C or S) or multi-element (C/S or C/H/S) analysis
- | Wide concentration range from 0.005 – 100 %
- | High weights possible for elemental analysis of heterogeneous samples
- | Optional autoloader with 36 or 130 positions (acid-resistant version available)




CARBON / HYDROGEN / SULFUR ANALYZER CHS-580A

## **OPERATION AND ANALYSIS PROCESS**

Operation of the CHS-580 A elemental analyzer is simple, fast and convenient. The samples are weighed and registered in the software as manual analysis or analysis via autoloader. Typical sample weights are between 50 and 500 mg. After starting the analysis, the sample is introduced into the furnace and the combustion gases CO<sub>2</sub> and SO<sub>2</sub> are measured in up to four infrared measuring cells. The results of the elemental analysis can be exported via LIMS, report or text file.



*Logging the sample into the software and weighing*

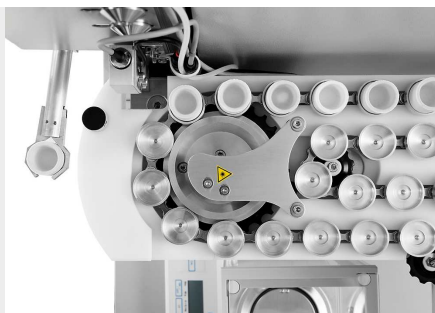

*Placing sample in autoloader and analysis*



*Measurement results and export*

## CARBON / SULFUR ANALYZER CHS-580A

### CONFIGURATIONS AND OPTIONS

ELTRA CHS-580 A analyzers can be configured to meet a wide range of customer requirements. C, H, S and C/ H/S configurations are available with up to 2 infrared cells for each element to cover a wide working range. Robust measuring cells for the reliable determination of sulfur in very high concentrations are also available.

The CHS-580 A series is perfectly equipped for TOC analysis according to DIN EN 15936 with optional acid-resistant autosamplers, as well as the optional TIC module.

An optional autoloader with 130 crucible positions is available for analyses with high sample throughput.



*Analyzer with optional TIC module*



*Autoloader with 130 crucible positions*



*Autoloader with acid-resistant plastic crucible holders*

CARBON / SULFUR ANALYZER CHS-580A

## TYPICAL SAMPLE MATERIALS

The CHS-580 A elemental analyzer is perfectly suited for measuring carbon, hydrogen and sulfur concentrations in matrices such as ashes, building materials, coal, coke, gypsum, limestone, oil, plant materials, rubber, sand, soil, soot, tobacco, waste, ...



*coal*



*soil*



*tobacco*

CARBON / SULFUR ANALYZER CHS-580A

## FUNCTIONAL PRINCIPLE

In the CHS-580A analyzers the sample is burnt in an oxygen atmosphere at temperatures of up to 1,550 °C. The furnace temperature can be freely selected in steps of 1 °C. The combustion gasses (CO<sub>2</sub>, H<sub>2</sub>O, SO<sub>2</sub>) coming from the furnace pass through a dust filter. After the water vapor is chemically absorbed, the dried CO<sub>2</sub> and SO<sub>2</sub> gas is detected in the additional infrared cells. Depending on the configuration, it is possible to combine up to four infrared cells with different sensitivities.

CARBON / SULFUR ANALYZER CHS-580A

## TECHNICAL DATA

|                               |                                                                                                                                             |
|-------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------|
| <b>Measured elements</b>      | carbon, hydrogen, sulfur                                                                                                                    |
| <b>Samples</b>                | organic                                                                                                                                     |
| <b>Furnace alignment</b>      | vertical                                                                                                                                    |
| <b>Sample carrier</b>         | ceramic crucibles                                                                                                                           |
| <b>Field of application</b>   | agriculture, biology, chemistry / plastics, coal / power plant, construction materials, environment / recycling, medicine / pharmaceuticals |
| <b>Furnace</b>                | resistance furnace (ceramic tube), adjustable up to 1550 °C (steps of 1 °C)                                                                 |
| <b>Detection method</b>       | solid state infrared absorption                                                                                                             |
| <b>Number of IR cells</b>     | 1 - 3                                                                                                                                       |
| <b>Material of IR path</b>    | gold                                                                                                                                        |
| <b>Typical analysis time</b>  | 60 - 180 s                                                                                                                                  |
| <b>Chemicals required</b>     | magnesium perchlorate, sodium hydroxide                                                                                                     |
| <b>Gas required</b>           | oxygen 99.5 % pure (2 - 4 bar / 30 - 60 psi)<br>compressed air (4 - 6 bar / 60 - 90 psi)                                                    |
| <b>Power requirements</b>     | 230 V, 50/60 Hz<br>max. heat up current 2000 W                                                                                              |
| <b>Dimensions (W x H x D)</b> | 55 x 100 x 60 cm<br>85 x 35 x 45 cm (Autoloader for 130 crucibles)                                                                          |
| <b>Weight</b>                 | ~ 90 kg (analyzer)<br>~ 15 kg (Autoloader for 130 crucibles)                                                                                |
| <b>Required equipment</b>     | PC, monitor, balance (resolution 0.0001g)                                                                                                   |
| <b>Required accessories</b>   | Autoloader (for 36 or 130 crucibles)                                                                                                        |
| <b>Optional accessories</b>   | TIC module, voltage stabilizer 5 KVA                                                                                                        |

[www.eltra.com/chs580A](http://www.eltra.com/chs580A)

## ORDER DATA

### ELTRA CHS-580A

(Please order PC, monitor, balance, Autoloader and consumables (starter-kit, anhydrous, sodium hydroxide) separately)

#### Measuring ranges at 500 mg sample weight

2)

|            |                                                                                    |          |                       |                                               |
|------------|------------------------------------------------------------------------------------|----------|-----------------------|-----------------------------------------------|
| 88100-4028 |   | CHS-580A | 1xH                   | 0.01 – 15% H                                  |
| 88100-4026 |   | CHS-580A | 1xC<br>+ 1xH          | 0.01 – 100% C<br>0.01 – 15% H                 |
| 88100-4027 |   | CHS-580A | 1xS<br>+ 1xH          | 0.005 – 2% S<br>0.01 – 15% H                  |
| 88100-4024 |   | CHS-580A | 2xC<br>+ 1xH          | 0.001 – 5% C   5% – 100% C<br>0.01% – 15% H   |
| 88100-4023 |  | CHS-580A | 1xC<br>+ 1xH<br>+ 1xS | 0.01 – 100% C<br>0.01 – 15% H<br>0.005 – 2% S |

Further measuring range combinations on request

### REQUIRED ACCESSORIES

#### PC, MONITOR, BALANCE

|            |                                                                                                                |
|------------|----------------------------------------------------------------------------------------------------------------|
| 71015-1000 | Computer with Intel Core i5-8400 Processor, 256 GB SSD; 8 GB RAM; Windows 10 operating system; keyboard; mouse |
| 88400-0584 | Monitor, TFT (23.8")                                                                                           |
| 88400-0645 | Balance (resolution 0.0001 g)                                                                                  |
| 14535      | Autoloader for 36 crucibles                                                                                    |
| 14535-1001 | Autoloader for 36 crucibles, acid resistant                                                                    |
| 14533      | Autoloader for 130 crucibles                                                                                   |
| 14533-1001 | Autoloader for 130 crucibles, acid resistant                                                                   |

#### REQUIRED CONSUMABLES / CHEMICALS FOR FIRST OPERATIONS

|            |                                                                                         |
|------------|-----------------------------------------------------------------------------------------|
| 88500-0006 | Starter-kit for 500 analyses CHS 580A (500 crucibles, 50 g glass wool, 100 g Combsolid) |
|------------|-----------------------------------------------------------------------------------------|

|       |                                                                                   |                                            |
|-------|-----------------------------------------------------------------------------------|--------------------------------------------|
| 90200 |  | Anhydron (magnesium perchlorate), 454 g 1) |
| 90210 |  | Sodium hydroxide, 500 g 1)                 |

## FURTHER OPTIONS AND CONSUMABLES

### ACCESSORIES (HARDWARE)

|            |                                                                                   |                  |
|------------|-----------------------------------------------------------------------------------|------------------|
| 88200-3800 |  | TIC-Module       |
| 72070      |                                                                                   | Oxygen regulator |
| 88400-0610 |                                                                                   | Barcode scanner  |
| 88600-0015 |  | Crucible holder  |

### CHEMICALS (FILLINGS FOR GLASS AND QUARTZ TUBES)

|            |                                                                                     |                                            |
|------------|-------------------------------------------------------------------------------------|--------------------------------------------|
| 90200      |  | Anhydron (magnesium perchlorate), 454 g 1) |
| 90210      |  | Sodium hydroxide, 500 g 1)                 |
| 90331      |  | Glass wool, 454 g                          |
| 90332      |  | Glass wool, 50 g                           |
| 92610      |  | Tube of high vacuum grease, 35 g           |
| 88600-0008 |  | Combsolid, 100 g                           |
| 91000-1005 |  | Copper, flakes, 25 g                       |
| 90840      |  | Quartz sand, 100 g                         |
| 88400-0508 |  | Steel wool                                 |

## CRUCIBLES

|            |                                                                                                                             |                                                              |
|------------|-----------------------------------------------------------------------------------------------------------------------------|--------------------------------------------------------------|
| 90149      |  | Ceramic crucibles, premium, Ø 1", foil-wrapped, 1,000 pieces |
| 90148      |        | Ceramic crucibles, premium, Ø 1", bagged, 1,000 pieces       |
| 88400-0176 |                                                                                                                             | Ceramic filtering crucibles, 100 pieces                      |
| 88400-0142 |                                                                                                                             | Filtering crucibles, 100 pieces, small version               |

## REAGENT TUBES

|       |                                                                                     |                                        |
|-------|-------------------------------------------------------------------------------------|----------------------------------------|
| 09090 |    | Reagent tubes 32 x 280 mm, 1 piece     |
| 09092 |    | Reagent tube XL size, 54 x 36 x 280 mm |
| 09093 |  | Reagent tube XXL size, length 280 mm   |

## TOOLS FOR OPERATION: SPATULAS, TWEEZERS, TONGS AND OTHER

|            |                                                                                     |                                                              |
|------------|-------------------------------------------------------------------------------------|--------------------------------------------------------------|
| 88400-0229 |  | Tweezers (160 mm), curved, 1 piece                           |
| 88400-0472 |  | Tweezers (145 mm), straight, 1 piece                         |
| 88400-0475 |  | Set with 6 spatula and 1 tweezers                            |
| 88400-0477 |  | Weighing boat, 1 piece, for weighing and usage of granulates |
| 90145      |  | Tongs for ceramic crucibles and boats, 220 mm, 1 piece       |
| 90146      |  | Steel pot for preheated crucibles and used crucibles         |

## TOOLS FOR MAINTENANCE

|            |                         |
|------------|-------------------------|
| 35400-8000 | Maintenance kit CS-580A |
|------------|-------------------------|

|            |                                                                                   |                                                                              |
|------------|-----------------------------------------------------------------------------------|------------------------------------------------------------------------------|
| 71010      |  | Brush, 16 mm, 1 piece, for cleaning balance from dust                        |
| 88400-0473 |  | Powder funnel (plastics), 1 piece, for easy filling of chemical tubes        |
| 88400-0490 |  | Rubber plug 29 x 35 x 30 mm, 1 piece, for sealing big glass tubes like 09090 |
| 88600-0026 |                                                                                   | Anhydrone filter tube                                                        |

## CALIBRATION MATERIALS

**Calibration materials may show slight variations depending on the current lot.**

**To see the current certification please visit [www.ELTRA.com](http://www.ELTRA.com).**

### COAL: CERTIFIED SULFUR CONTENT

|            |                                                                                     |                           |
|------------|-------------------------------------------------------------------------------------|---------------------------|
| 92511-3005 |                                                                                     | Coal, 50 g, < 0.1 % S     |
| 92511-3010 |   | Coal, 50 g, 0.1 – 0.5 % S |
| 92511-3020 |  | Coal, 50 g, 0.5 – 1.0 % S |
| 92511-3030 |  | Coal, 50 g, 1.0 – 1.5 % S |
| 92511-3040 |  | Coal, 50 g, 1.5 – 2.0 % S |
| 92511-3050 |  | Coal, 50 g, 2.0 – 3.0 % S |
| 92511-3060 |  | Coal, 50 g, 3.0 – 4.0 % S |
| 92511-3070 |  | Coal, 50 g, 4.0 – 5.0 % S |
| 92511-3080 |  | Coal, 50 g, >5.0 % S      |

### COAL, PREMIUM, C/H/N/S, ASH, VOLATILE CONTENT, CERTIFIED

|            |                                |
|------------|--------------------------------|
| 92550-3010 | Coal, premium, 50 g, < 1.0 % S |
|------------|--------------------------------|

|            |                                                                                   |                                    |
|------------|-----------------------------------------------------------------------------------|------------------------------------|
| 92550-3020 |  | Coal, premium, 50 g, ~ 1 % S       |
| 92550-3040 |  | Coal, premium, 50 g, 1.0 – 3.0 % S |
| 92550-3060 |  | Coal, premium, 50 g, > 3.0 % S     |

#### COKE, PREMIUM, C/H/N/S, ASH, VOLATILE CONTENT, CERTIFIED

|            |                                                                                   |                     |
|------------|-----------------------------------------------------------------------------------|---------------------|
| 92560-3010 |  | Coke, premium, 50 g |
|------------|-----------------------------------------------------------------------------------|---------------------|

#### PET COKE, PREMIUM, C/H/N/S, ASH, VOLATILE CONTENT CERTIFIED

|            |                                  |
|------------|----------------------------------|
| 92570-3020 | Pet coke, premium, 50 g, ~ 1 % S |
| 92570-3040 | Pet coke, premium, 50 g, ~ 1 % S |

#### OTHER

|            |                                                                                     |                           |
|------------|-------------------------------------------------------------------------------------|---------------------------|
| 90812-3001 |  | Limestone, 25 g, 0.04 % S |
|------------|-------------------------------------------------------------------------------------|---------------------------|

|            |                                                                                     |                          |
|------------|-------------------------------------------------------------------------------------|--------------------------|
| 90812-3002 |  | Limestone, 25 g, 0.4 % S |
|------------|-------------------------------------------------------------------------------------|--------------------------|

|            |                          |
|------------|--------------------------|
| 90812-3003 | Limestone, 25 g, < 5 % C |
|------------|--------------------------|

|            |                             |
|------------|-----------------------------|
| 90812-3004 | Limestone, 25 g, 5 – 10 % C |
|------------|-----------------------------|

|            |                        |
|------------|------------------------|
| 90817-3001 | Soil, 25 g, > 2 % C, S |
|------------|------------------------|

|            |                        |
|------------|------------------------|
| 90817-3002 | Soil, 25 g, < 1 % C, S |
|------------|------------------------|

|            |                        |
|------------|------------------------|
| 90817-3003 | Soil, 25 g, > 2 % C, S |
|------------|------------------------|

|            |                        |
|------------|------------------------|
| 90817-3004 | Soil, 25 g, > 2 % C, S |
|------------|------------------------|

|            |                                                                                     |                     |
|------------|-------------------------------------------------------------------------------------|---------------------|
| 91900-1001 |  | Ore, 30 g, ~1.4 % S |
|------------|-------------------------------------------------------------------------------------|---------------------|

|            |                                                                                     |                     |
|------------|-------------------------------------------------------------------------------------|---------------------|
| 91900-1002 |  | Ore, 30 g, ~4.2 % S |
|------------|-------------------------------------------------------------------------------------|---------------------|

|            |                   |
|------------|-------------------|
| 91900-1003 | Ore, 30 g, ~3 % S |
|------------|-------------------|

#### PRIMARY SUBSTANCES FOR CALIBRATION

90710-3010



EDTA, 50 g

90710-3030

Sucrose, 50 g

90810



Calcium carbonate, 100 g

90821



Barium sulphate, 50 g

90824



Sulfanilic acid, 50 g

91900-2001

Zinc sulfide, 50 g, 32 % S

**Please note: Every analyzer requires PC, monitor, balance and some consumables (crucibles, chemicals) which have to be ordered separately**